ARREDAMENTI  
PER LABORATORI

**MA95**  
**EVO** 

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**TALASSI**

## TECHNICAL FURNITURE FOR LABORATORIES

Talassi designs and manufactures technical furniture for laboratories since 1969. Today it represents an important landmark in the world of scientific laboratories: more than forty years of experience serving the national and international scientific community with a wide range of products. The dimension and organization of the Company allow to keep flexible the processes of designing and manufacturing adapting the production to the customer needs. Reduced construction time and customer tailored production differentiate Talassi from the competitors, more accustomed to massive and standardized productions. Our technical furniture complies with the specific needs of chemical-pharmaceutical, microbiological and instrumental analysis laboratories. Our customers are steel industries, food industries, manufacturing industries, chemical and pharmaceutical industries, research laboratories, hospital, public and private laboratories, technical schools and governmental scientific institutes of control. We strongly believe that to design a practical and comfortable laboratory needs the full understanding of the procedures and methods of the end user. The creation of a project requires a deep harmony with our client enabling him to share with us his own experience. A safe laboratory is a balance between the guidelines and a proportioned distribution of the furniture. Every year we invest in research and development to create new solutions and integrate new technologies into our products. We give great importance to the design of new products and to the research of new materials able to limit the environment impact favouring recyclable materials use : metal, glass and wood. Talassi is certified UNI EN ISO 9001 and UNI EN ISO 14001 The furnishings is certified EN 13150 and EN 14175. Talassi is registered to the main platforms for the electronic market of the public administration: SINTEL and MEPA.

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## SAFETY CHEMICAL HOOD

*The safety in the laboratory depends on combined factors among which a crucial role is played by the knowledge and the use of collective protection systems (CPD). The fume hood ensures the operator protection during the processing of potentially hazardous materials, is defined therefore as a Collective Protection Device. The choice of the chemical hood so it is of paramount importance in the context of a laboratory, you must assess all the risks faced by the operators, the frequency of use of hazardous substances, the level of security required, the lay-out of the laboratory, the ergonomics and last but not least the energy saving. A careful study of these factors along 45 years of experience in the preparation of scientific laboratories in close contact with end users has led us to the creation of the new chemical fume hood MA95EVO.*

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CERTIFICATION  
CERTIFICACIÓN

**PANNELLO DI CONTROLLO  
MA95 FLOW**

MA95 FLOW CONTROL PANEL  
PANEL DE CONTROL MA95 FLOW

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BOARD WITH SOCKETS  
PANEL CON ENCHUFES

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## THE EVOLUTION OF THE FUME HOOD

MA95 EVO is the evolution in construction and functional philosophy of the fume hoods, it represents the state of the art in terms of performance and safety. The traditional shape includes new contents, with particular attention to the study of aerodynamic flows and electronic controls that assure increased aerodynamic performance and an unprecedented energy saving. The chemical hood MA95 EVO is the synthesis of the control of the flow of air for the total protection of the user in any operating condition. The careful study of aerodynamics guarantees to the hood a modern design and high-performance in operational functionality, endorsed by the report certification according to the standards EN14175. The particular internal aerodynamics conformation has allowed to achieve excellent performance of containment of sulphur hexafluoride SF<sub>6</sub> with values <0.1 ppm even at reduced air flow (as evidenced by the test report according to EN 14175-3-6) allowing an appreciable saving in energy consumption. The ergonomic position of the commands of the technical services allows the operator easy manipulation in safety conditions. The MA95 EVO is also available in walk-in configuration and with services commands installed along the sides of the fume hood.

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## **EVO MA95: STANDARD EQUIPMENT**

- Supporting structure in welded steel painted with epoxy powder in heated tunnel
- External sides in steel sheet, painted with epoxy powder in heated tunnel
- Sash sliding vertically
- Protective safety glass sliding horizontally
- Electronic control MA95FLOW with display and visual and sonorous alarm
- Multilingual software
- Door locking device at 50 cm opening
- SAFETY-BLOCK device in case of ropes breaking
- Fume hood plenum with vents, removable for inspection and maintenance
- Worktop in monolithic GRES with retaining edges, sink and syphon
- Internal channel to carry moisture to the worktop
- Equipped to connect stainless steel framework
- Saving energy Led lamp IP65 for indoor lighting > 400 lux
- Lines services: water, gas
- IP55 220V electrical outlets protected by circuit breaker
- Undertop cabinet with hinged doors

## **ACCESSORIES AND / OR CUSTOMIZATIONS**

- Box with activated charcoal filters
- Box with combined charcoal and absolute filters
- Cabinets for acids and bases
- Cabinets for flammables
- Disposal containers for dangerous liquid
- Internal coatings (high-resistance polypropylene, PVC, stainless steel, stoneware)
- Stainless steel frameworks for glassware
- Technical gases taps
- Fume hood models with services control on the sides and Walk-in configuration
- Worktops in AISI 316 stainless steel, polypropylene or Corian

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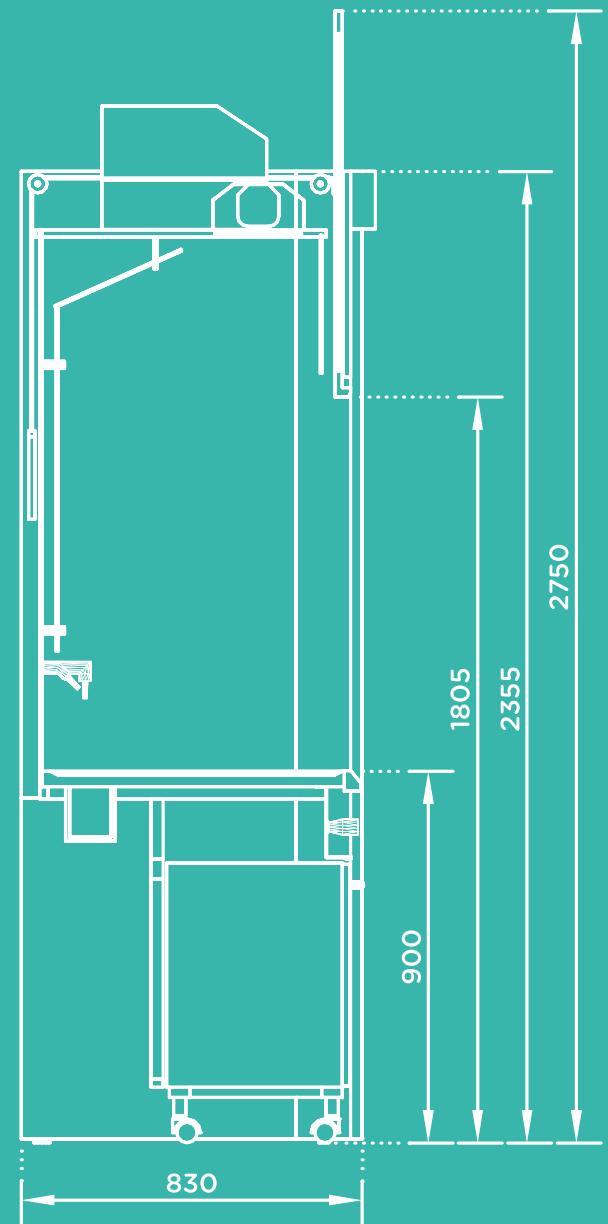
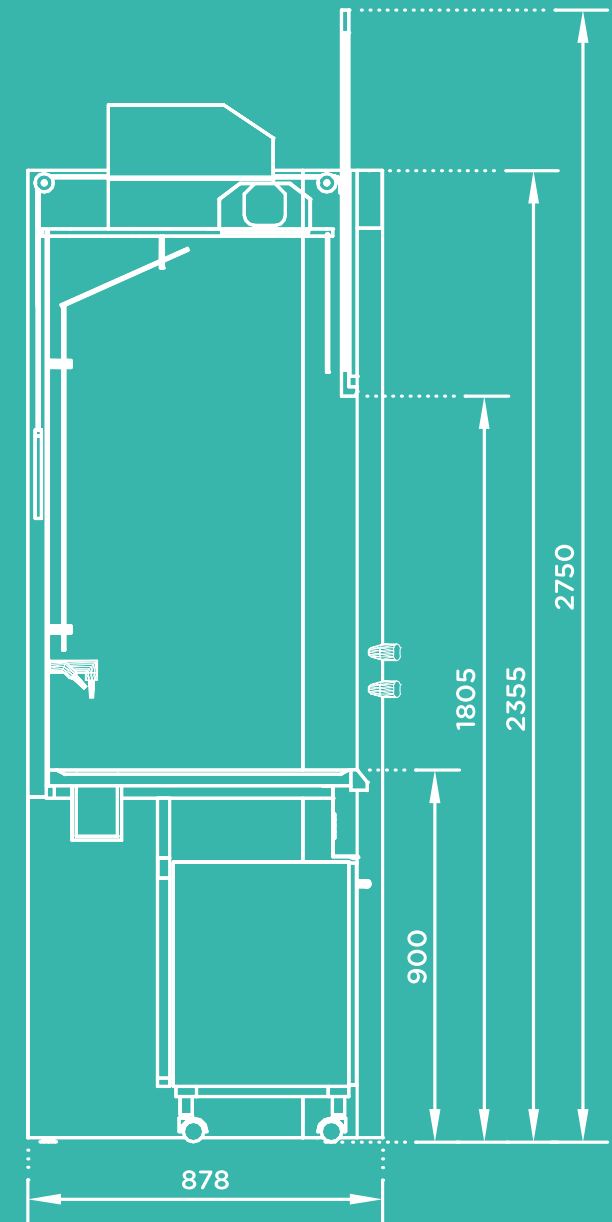
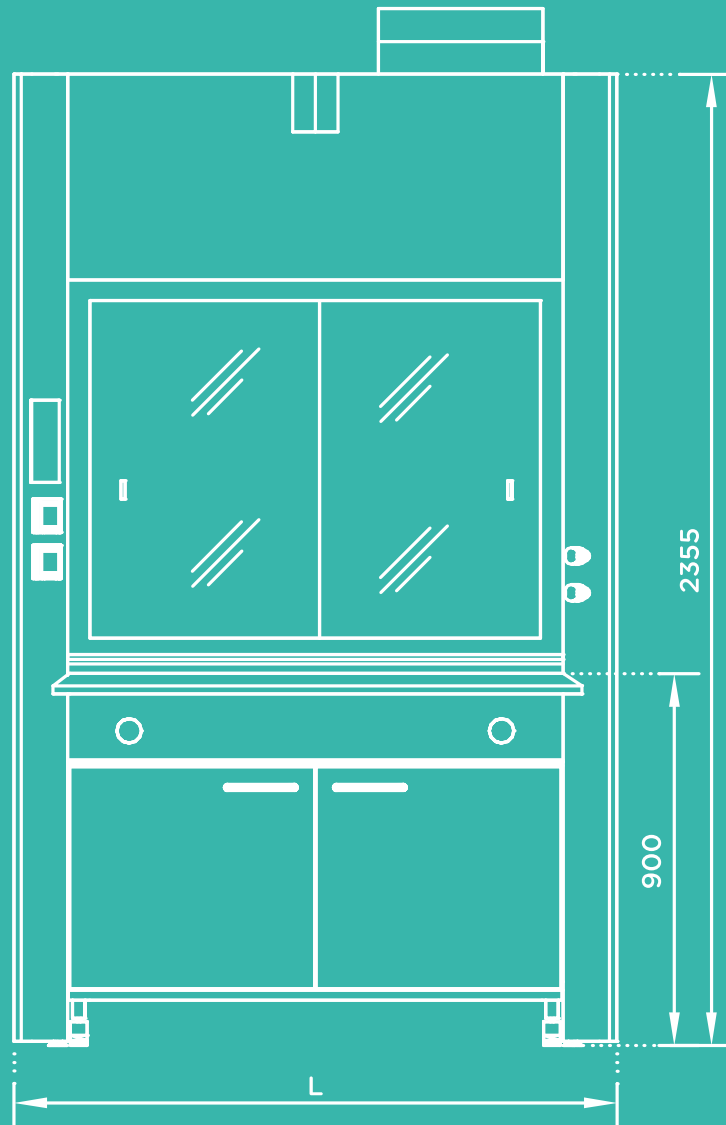
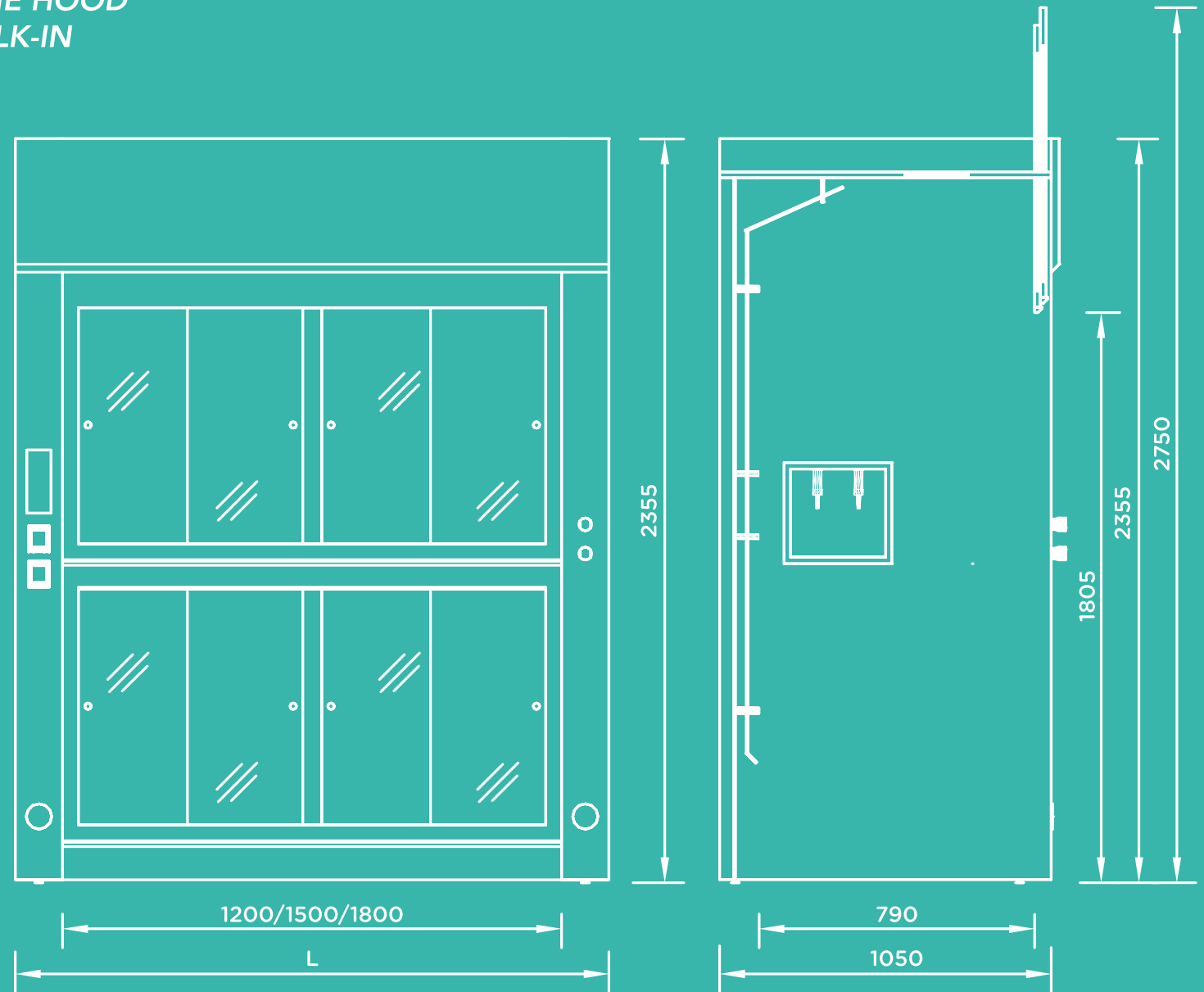
### MODELLO CON SERVIZI SUI FIANCHI / MODEL WITH SERVICES CONTROL ON THE SIDES

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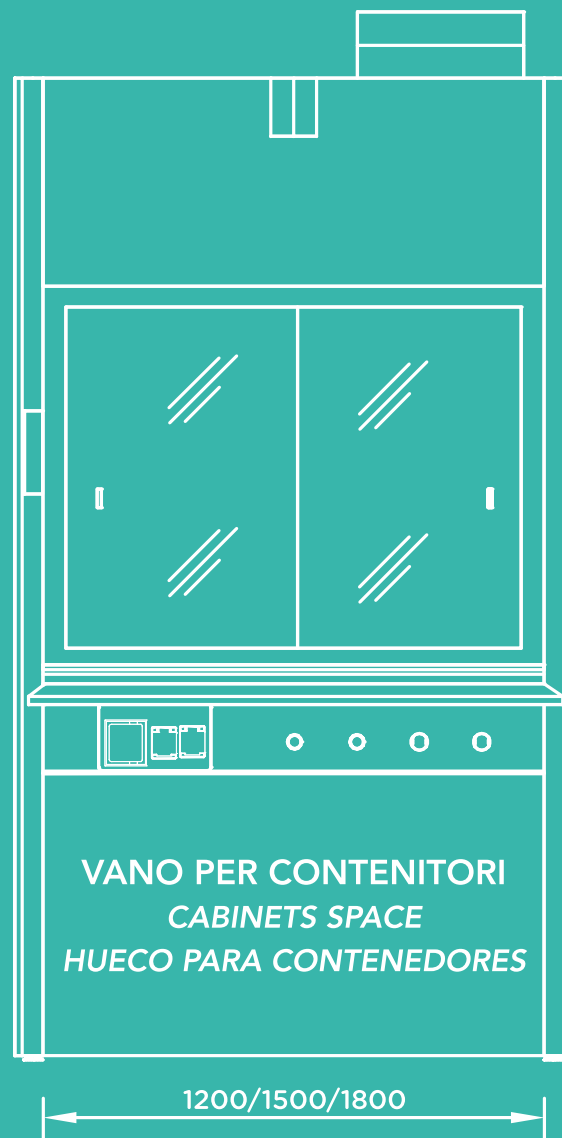
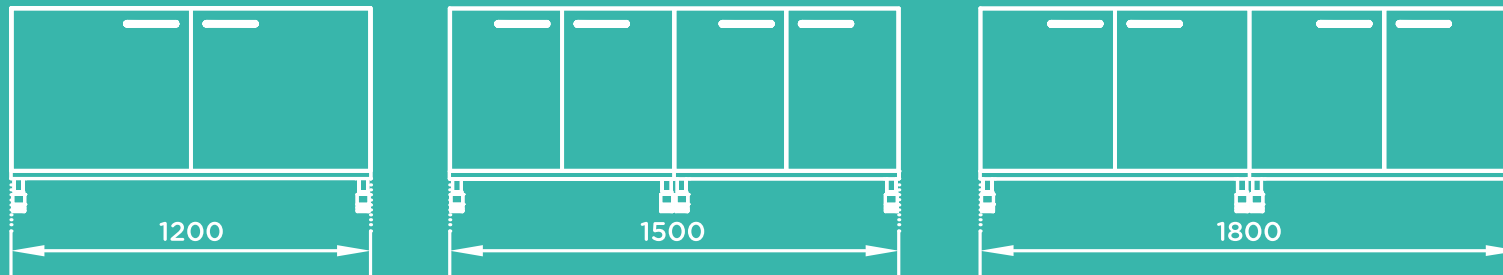
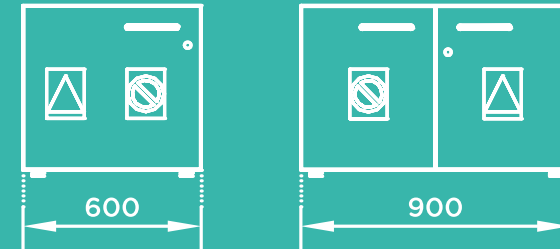
*CONTENEDORES PARA INFLAMABLES*


## CONTENITORI ASPIRATI FILTRATI

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*ASPIRATED AND FILTERED CABINETS FOR ACIDS AND BASES*

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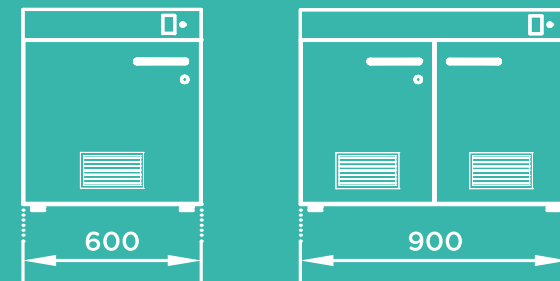### MA95 FLOW

*The electronic control MA95FLOW, with screen display, is an intuitive interface that allows to monitor in real time all the parameters of operation of the fume hood assuring an high safety standard. Multilingual programming, operation detector, visual and audible alarm in case of insufficient suction or in case the sash exceed 50 cm height opening, control of starting, internal lighting control, hot-wire probe management system, maintenance service management . Equipped to control inverters, motor valves, motorized sash. The system can be easily implemented even in fume hood already installed.*

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TALASSI  
BIO-EDGE EOP 4  
1102200  
1100  
Trade by  
AUTO & SAFE  
MAX  
MAX 90 FLOW

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## AUTOMATIC FLOW DRIVE

*The fume hood MA95 EVO has been equipped with a valuable accessory which automatically adjust the airflow (VAV). When the fume hood is operating in automatic mode the control system MA95 Flow monitors and adjusts the air intake flow in relation to the opening of the sash. This regulation, known as "variable flow", ensures the minimum speed able to assure the containment of toxic substances near the sash. The traditional fume hoods do not have the ability to adjust the volume of air extracted from the room where they are installed so the energy consumption is constant and therefore can not be decreased when the sash is lowered. This means waste of electricity and, especially, consumption above the real needs of conditioned air. The fume hood MA95 EVO equipped with the VAV system is able to maintain the necessary level of security, reducing at the same time consumption of, heated or cooled, air and of electrical energy for the fan engine.*

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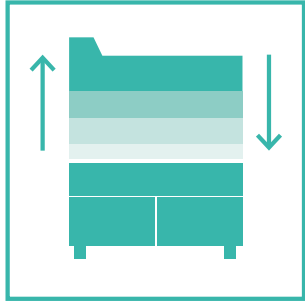
## AUTOMATIC SASH CLOSING

*The self-closing sash is more and more required by laboratory technicians. The hood MA95 EVO can be equipped with a device that, after a scheduled time, automatically close the sash. The sash opening is the focal point of the protection of the hood, but also the element that, if closed, lowers the consumption of air drawn from the laboratory. In fact if the automatic flow drive (VAV) drastically decreases the air flow rate when the sash is closed, the automation ease the systematic closing of the same. The automation of the sash grants (in the absence of the operator) the full shut down. When the sash is closed, the vacuum flow from the fume hood, thanks to the VAV, is reduced up to 75%. The coupling of the VAV system with the automatic sash closing system increases the energy saving but above all, preventing accidents, the laboratory safety.*

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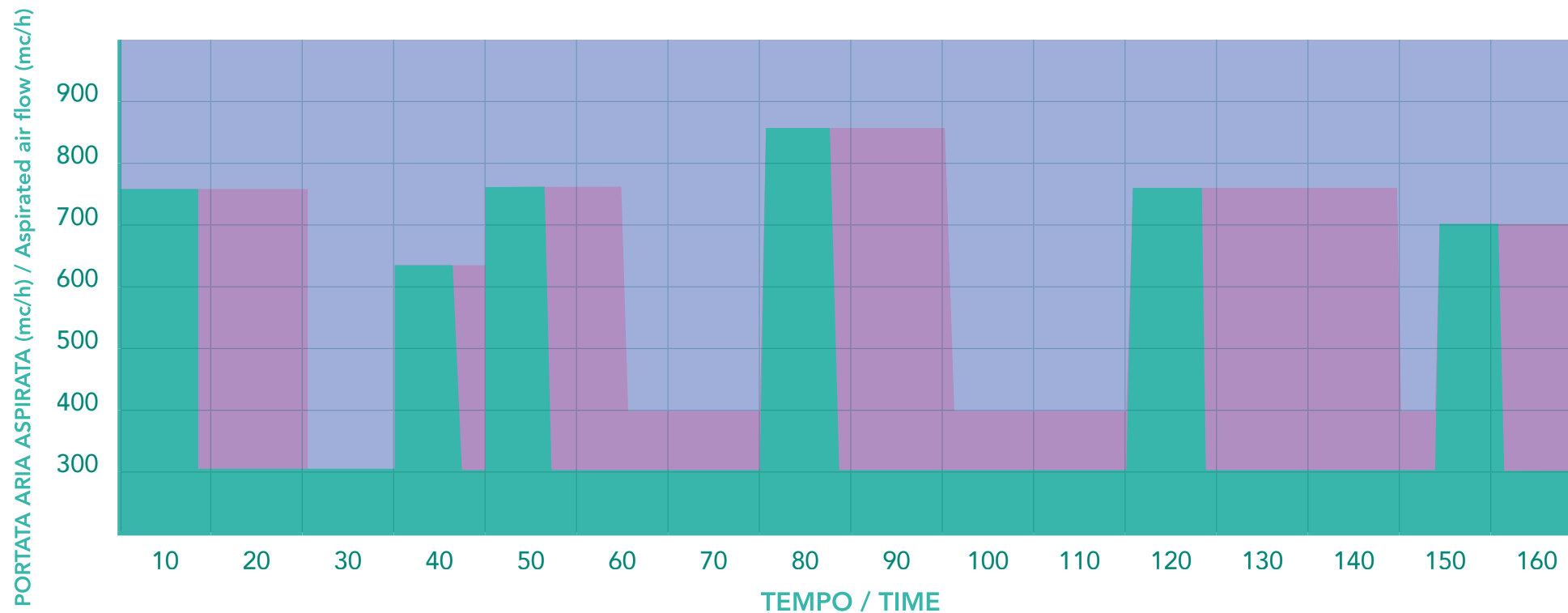
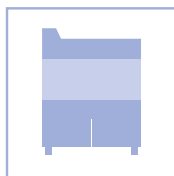
AIR EXTRACTED FROM  
THE FUME HOOD WITH  
CONSTANT FLOW

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AIR EXTRACTED FROM  
CHEMICAL HOOD WITH  
VARIABLE FLOW

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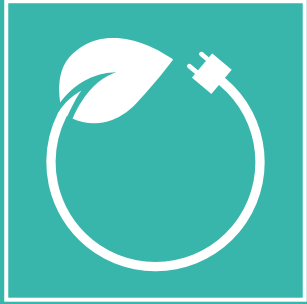
AIR EXTRACTED FROM CHEMICAL  
HOOD WITH VARIABLE FLOW AND  
AUTOMATIC CLOSING SASH

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### **Comparison of energy consumption**

*In the diagram examples are indicated the values of the average consumption of a fume hood, 120 cm wide, in the three possible configurations in relation to the working time.*

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## CHARACTERISTICS AND CERTIFICATIONS

- *Quality system certification according to UNI EN ISO 9001*
- *Environmental system certification according to UNI EN ISO 14001*
- *Product certification of chemical hoods UNI EN 14175-1-2-3-6*
- *Product certification of benches UNI EN 13150*
- *Product certification of laboratory cabinets UNI EN 14727*
- *Product certification of cabinets for flammable products UNI EN 14470-1*
- *Product certification for electrical safety in laboratory UNI EN 61010-1*

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### MAINTENANCE SERVICE

*In addition to the design, production and installation, Talassi offers a qualified maintenance service of the fume hoods and tests on site according to UNI EN 14175-4.*

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